



Valiteq Media Fill Test Kit

The AAOA has worked to develop resources and tools to help you meet the compliance criteria outlined in the pending chapter.

Last Fall, we released our [USP General Chapter <797> Sterile Allergen Extract Compounding Online Module](#) to help you and your practice meet the compounding personnel criteria. USP <797> requires personnel to demonstrate proficiency in knowledge of principles and skills for sterile compounding every 12 months and to document training and competency. Partnering with Area9 and its adaptive learning platform, the AAOA's online module is a great tool to help assure you can meet and document the proficiency defined in USP <797>. Because it is based on adaptive learning, module participants will succeed through the learning process. So, it is a great tool to help learn about sterile compounding as well as document proficiency.

AAOA has also partnered with [Valiteq](#) to offer kits (AllerTEQ A1CDI) at a special AAOA member rate to meet the media fill and gloved fingertip testing requirements. AAOA members will get a 15% discount on the kits. The kits list for \$200 and AAOA members will pay \$170 plus shipping.

While the implementation dates for the new USP General Chapter <797> are still undefined, AAOA's [resources](#) are available for you and your staff as you position your practice to meet the new guidance.

Here are some step-by-step instructions on how to order your test kits and how to handle and store it once received.

To order your test kits, please fill out the attached form and use the code AAOAUSP2020. The form should be emailed to aaoa@valiteq.com.

Each kit, currently defined as AllerTEQ A1CDI, consists of the media fill test, the gloved fingertip test kit with 6 contact plates, 2-week incubation process, 1-week growth promotion, and a written report at 4 weeks.




Valiteq Media Fill Test Kit

Receiving and handling your test kits

Your shipment arrives by second day air and it contains perishable contents, so it's important to follow the instructions on how to unpack, store and return the test kit.

Notice - Do not discard any of the contents, until the test is completed and ready to be returned Valiteq for analysis.

1. Content sheet

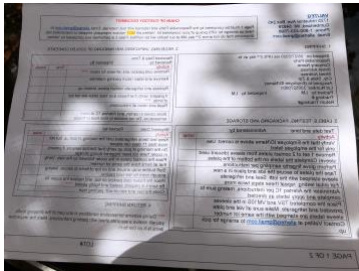

- Describes what is in the test
- Denotes employee doing the test
- Initial each activity

This form needs to be returned along with the test once the test is completed.

2. Unpacking Instructions



Includes unpacking instructions for normal weather conditions as well as for warm weather shipping & packaging (plates need to be refrigerated). Please read the 15 steps before unpacking and storing your test kit to ensure unpacking and storing is done correctly. Also, you'll find the return instructions.

3. Test kit and plates:




Your shipment contains two boxes. The smaller one is the Media fill test kit. The larger box labeled Perishable contains the 3 plates (follow instructions on how to correctly store).



The American Academy
of Otolaryngic Allergy

Valiteq Media Fill Test Kit

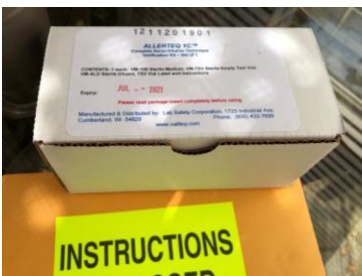

Media fill test kit (vials, instructions and shipping information)



3 Testing Plates – Need to be refrigerated

If you have any questions, please check out our Unboxing Video or call our office at 202.955.5010.